

SPOTCHECK™ & SPOTCHECK PLUS™ Non Instrument Tests

SPOTCHECK™-Glucose Test

SPOTCHECK™ detects Glucose, present in 85% of food products, without instrumentation. The self-contained reagent turns from clear to green when food residue is present in a surface sample. With SPOTCHECK™, you detect almost instantly if a surface is unclean, allowing immediate corrective action.



SPOTCHECK PLUS™-Lactose & Glucose Test

All the benefits and convenience of Glucose detection from SPOTCHECK™ with an added Lactose detection reagent. Perfect for use in the dairy industry.




Key Features

- Results in under 60 seconds
- Detection comparable to ATP systems
- All-in-One Device - No Instrumentation
- Allows for immediate corrective action
- Easy to interpret
- Affordable

Ideal for Processors & Handlers of:

- Baked Goods
- Fruits & Vegetables
- Beverages
- Dairy Products
- Confectionary Products
- Dressings & Sauces

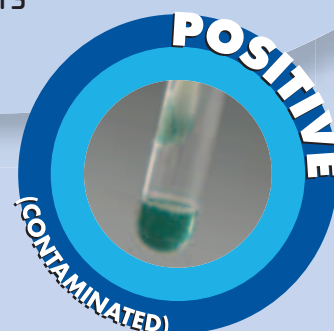

hygiena

Rapid Solutions for Food Safety

941 Avenida Acaso, Camarillo, CA 93012

tel: 805.388.8007 ext. 300

e-mail: info@hygienausa.com

web: www.hygienausa.com

Comparative Performance Sheet

Industrial Food Processing

Lowest Dilution Detected

SAMPLE TYPE	COMMERCIAL PROTEIN TEST	SPOTCHECK™ & SPOTCHECK PLUS™	ATP BIOLUMINESCENCE
Dairy Products	1/10	1/100	1/100 - 1/1000
Meat	1/10	1/10	1/100
Root Vegetables	Undetected	1/10	1/50
Salad Vegetables	Undetected	1/100	1/100
Fruits	Undetected	1/500	1/500
Bakery Products	Undetected	1/10	1/100
Confectionary	Not Tested	1/10 - 1/1000	1/10 - 1/1000

Food Service/Catering

SAMPLE TYPE	COMMERCIAL PROTEIN TEST	SPOTCHECK™ & SPOTCHECK PLUS™	ATP BIOLUMINESCENCE
Salad Vegetables			
Lettuce	Undetected	1/10 - 1/100	1/1000
Tomato	Undetected	1/10	1/10000
Pickle	Undetected	Undetected	1/100
Dressings			
Salad Dressing	1/10*	1/10 - 1/100	1/10
Ketchup	1/10*	1/1000	1/100
Beverages			
Soft Drinks	Undetected	1/1000	1/1000
Fruit Juices	Undetected	1/1000	1/10000
Milk Shake	1/10*	1/100	1/100
Bakery Products			
Bread	1/10	1/100	1/100
Pie	1/10*	1/100	1/100

Data presented equals the minimum dilution of the product type successfully detected by each method (at least three different food types in each category, and three replicates per test).

* Color change was extremely slight



With SPOTCHECK™ & SPOTCHECK PLUS™, you can detect all major food groups, unlike protein tests that are limited to high-protein foods such as meat.

Contact us for free samples.



hygiene

Rapid Solutions for Food Safety

941 Avenida Acaso, Camarillo, CA 93012
 tel: 805.388.8007 ext. 300
 e-mail: info@hygieneusa.com
 web: www.hygieneusa.com