


How to Change Your Alarm Settings (FridgeTag)

8.1. Initial menu (read and change settings)

To change the date format, the date, the time, the temperature measurement unit or the alarm settings or to read the preset alarm limits please proceed as follows:

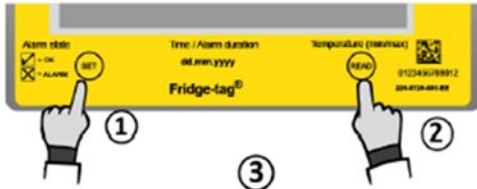

1. Press and hold SET ...
2. ... then press READ shortly ...
3. ... then release both buttons simultaneously.

SET DATE (internal sensor) is now displayed on the screen.

OUTSIDE (external sensor) is now displayed on the screen.

You entered the menu mode and may choose which entry to see or change.

You can access the following 4 menus:

OUTSIDE (external sensor): first screen, shows the temperature measured with the internal sensor of the Fridge-tag (normal ambient temperature).

Press READ once to get to SET DATE.

SET DATE (internal sensor): Configuration with internal sensor,
SET DATE is directly shown.

1. SET DATE: change date and/or time settings
2. READ CONF: read the alarm settings
3. CELS FAHR: change the temperature unit
4. SET CONF: change the alarm settings (only if enabled in factory setting)

Use the READ button to scroll through the menu.

Use the SET button to access the corresponding menu.

Select 4. SET CONF

For Fridge

- **0:05, 8.0** (set to alarm when warmer than 8°C for longer than 5 minutes)
- **0:10, 2.0** (set to alarm when colder than 2°C for longer than 10 minutes)

For Freezer

- **0:05**, (after entering 0:05 press set one time, press read until you see a negative symbol, then press set) **-15** (alarm when warmer than -15°C for longer than 5 minutes)
- **0:10**, (after entering 0:10 press set one time, press read until you see a negative symbol, then press set) **-25** (alarm when colder than -25°C for longer than 10 minutes)