

STEELCO LAB 610: QUICK-START GUIDE


STEP 1: PRE-WASH CHECK




Wear required PPE



Remove all labels and tape from glassware

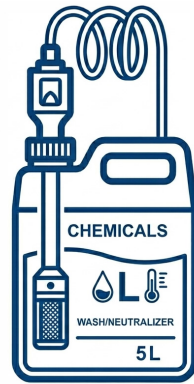

Confirm chemical levels.
Check for 'No/Low Chemical alarms'

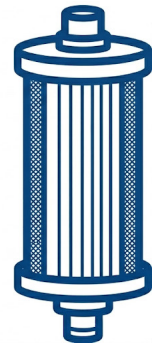


Clear debris from the filtration system

STEP 2: RACK SELECTION



Use for beakers, jars, misc. items

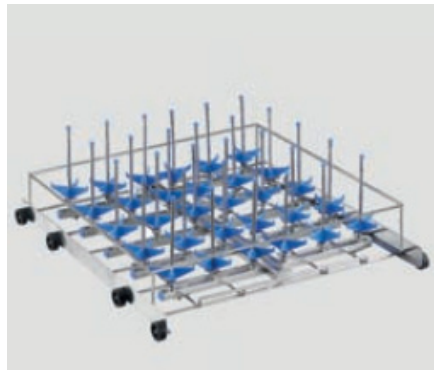

Use for narrow-neck flasks/cylinders



Use for pipettes
(point tips down)

STEP 3: LOADING & TROLLEY

Note:

Place glassware face down and ensure items are secured in place when loading


Trolley Instructions:

1. Align trolley with open washer door to dock
2. Engage safety latch to lock trolley to washer
3. Slide rack from trolley into the washer chamber till it is aligned with the water injectors
4. Disengage trolley and close door



STEP 4: START CYCLE

1. Turn knob to ON position
2. Press **P+** to select a program
3. Verify program on screen
4. Press **⏻** to start cycle

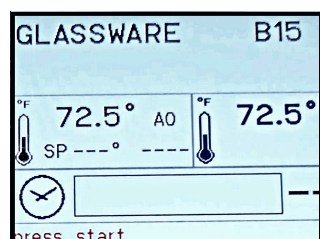

ALARM RESET

1. Hold **START** **⏻** + **STOP** **⏻** simultaneously for 5 seconds
2. Press **P2** once, then **P1** once to clear alarm