


Model AF Series Lab Ovens

OPERATING MANUAL

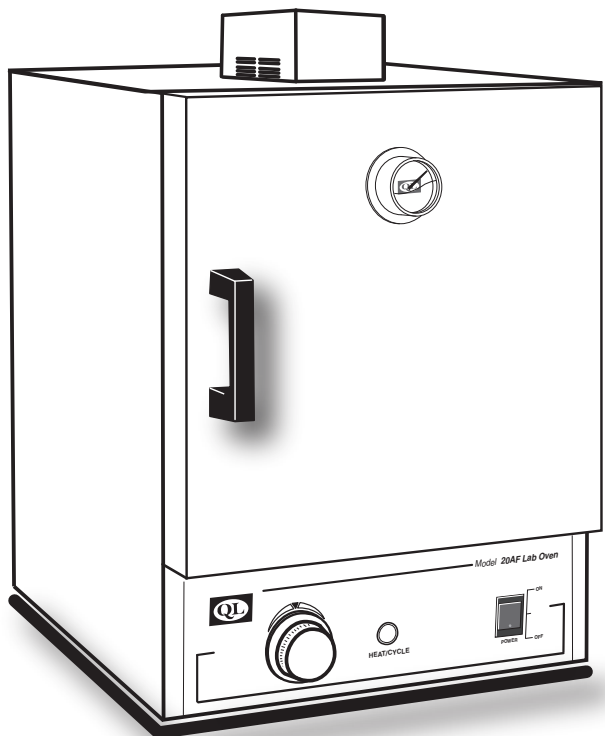

Carton Contents

- (1) AF Series Oven
- (2) Adjustable chrome wire shelf
- (4) Shelf brackets

Approvals:

Underwriter's UL / CUL, United States/
Canadian for laboratory equipment.

Compliance: UL Standard 61010-1
IEC 61010-1.

NOT FOR USE WITH FLAMMABLE LIQUIDS OR GASES



SPECIFICATIONS	MODEL 10AF	MODEL 20AF	MODEL 30AF	MODEL 40AF
INTERIOR DIMENSIONS				
INCHES W x H x D	12x8.25x10	13x11x13	18x14.13x12	18x19.63x14
(CM) W x H x D	30.5x21x25.4	33x28x33	45.7x35.5x30.5	45.7x50x35.5
EXTERIOR DIMENSIONS				
INCHES W x H x D	14x20.5x12.25	15x25x15.25	20x29x14.25	20x31.5x16.25
(CM) W x H x D	35.5x51x31	38x63.5x38.75	51x73.7x36.2	51x80x41.25
WEIGHT (lbs)				
SHIPPING	50	66	82	112
STAND ALONE	44	49	74	95
TEMPERATURE RANGE				
Ambient + 25°F to F / C	450°/232°	450°/232°	450°/232°	450°/232°
TEMPERATURE STABILITY				
Typically +/- F / C	6° / 3°	6° / 3°	6° / 3°	6° / 3°
STANDARD ELECTRICAL*				
VOLTS/AMPS	120/6.7	120/8.8	120/12.5	120/12.5
FREQUENCY/HZ	50/60	50/60	50/60	50/60
WATTS	800	1050	1500	1500
PLUG/NEMA	5-15P	5-15P	5-15P	5-15P

* Standard models voltage only, optional 220 voltage available. For power requirements, see electrical info on label located at the rear of the oven.

Common Unit Construction

Exterior: Powder-Coated Steel







Insulation: Fiberglass

Thermo-control: Bimetal/Hydraulic


Interior: Aluminized Steel

Motor: Sleeve Bearing, Thermally Protected


Heater: Resistive-Tubular Incoloy

-  **WARNING:** READ INSTRUCTIONS THOROUGHLY BEFORE OPERATING OVEN
-  **WARNING:** RISK OF ELECTRICAL SHOCK. DISCONNECT UNIT FROM POWER SOURCE BEFORE REMOVING COVER
-  **CAUTION:** EXTERIOR SURFACES MAY BECOME HOT DURING OPERATION
-  **WARNING:** DO NOT USE WITH FLAMMABLE LIQUIDS OR GASES
-  **WARNING:** NOT FOR USE WITH OPEN LIQUIDS
-  **WARNING:** CARE SHOULD BE TAKEN WHEN CARRYING/LIFTING UNIT INTO PLACE. LARGER UNITS ARE HEAVY AND MAY REQUIRE TWO PEOPLE TO TRANSPORT OR INSTALL.

Safety Precautions  Read Operating Instructions Thoroughly Prior to Operation

 The AF series Lab Ovens are not designed for use with any flammable solvents or gases or for materials that may contain flammable solvents or gases, or in close proximity to any flammable solvents, gases or materials. Operate oven in accordance with all state, local, and federal laws that may govern its use for your specific application. Use only a grounded outlet that is rated for your model's electrical requirement. Oven exterior walls and doors may become hot to the touch when operating at higher set temperatures. Do not leave the oven unattended during operation, especially when processing materials that have flash point temperatures lower than the oven model's maximum operating range. Do not modify the oven or factory control settings to operate the oven above the stated maximum operating temperature. Conduct periodic maintenance as required.

Set-up & Installation

Set-up unit in a clean and unobstructed operating location. Keep a minimum of 4 inches of airspace around the unit and a minimum of 18 inches above the unit. The 2 exhaust ports at the top of the unit will expel a small amount of heated air through forced convection. This "fixed" amount of exhausted air gently removes any fumes or moisture from the chamber during processing. This prevents condensation within the chamber and prolongs the chamber life over time.  Do not place anything on top of the motor cover. This will severely overheat the motor and cause risk of fire. **(FIG 1)**

Important: The exhaust ports should NOT be used as chamber access for temperature-measuring probes. Insertion of any such probe or device may damage or unbalance the circulating fan blade at the top of the oven chamber.

Plug oven into a grounded outlet rated for your unit's labeled voltage. A separate circuit or breaker should be dedicated for the oven.

Install adjustable shelves as required, by placing the ends of the wire shelf bracket into the corresponding holes located on the inner sides of the oven at the desired height. Push ends of the bracket into the holes until the first bends in the bracket are against the wall, then rotate the bracket down. Place the shelf on the brackets. **(FIG 2)**

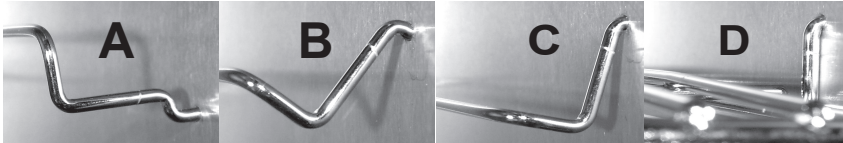


FIG. 2

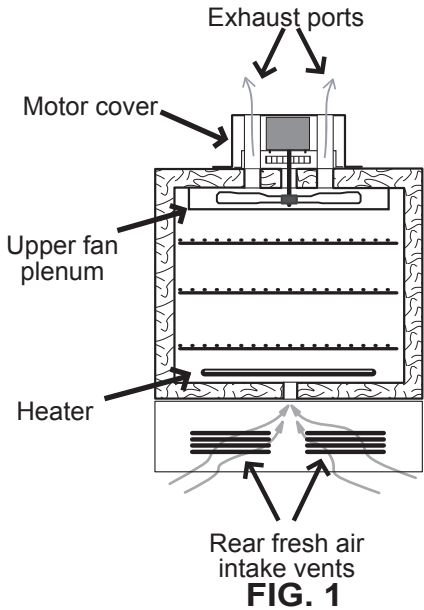


FIG. 1

General Operation

All models are ready for your immediate use. Turn the ON/OFF power switch to the ON position. The fan will begin to circulate air and the oven will begin to heat toward a temperature setting.

To set the temperature for the 10AF series model: Turn the thermostat control knob clockwise to around the number '8'. The heat-cycle light will come on indicating heater activity. Monitor the reading of the door thermometer until it reaches the desired temperature. Next, rotate the control knob very slowly counter-clockwise until the heat cycle light turns off. The thermostat will then cycle automatically, maintaining the set temperature. Turn oven off with the ON/OFF power switch, saving the temperature setting for the next use.

For 20AF, 30AF, and 40AF models: These models come equipped with a dual-range "quick-set" knob dial. This feature makes it easier to hit a target temperature range from a cold start, whether using the Fahrenheit or Centigrade scale. Turn the thermostat knob to the desired temperature indicated on the knob (F), or the dial (C).



FIG. 3

To set a temperature in degrees Fahrenheit, turn the knob in a clockwise direction until the desired number on the knob is directly under the white triangle with the "°F" marking. **(FIG 3)**

General Operation (continued)

To set a temperature in degrees Centigrade, turn the knob in a clockwise direction until the white triangle with the "°C" marking is pointing toward the desired temperature printed on the face of the control panel. (FIG 4)

Rotate the thermostat dial to the desired temperature. The heat-cycle light will illuminate until the set temperature is reached. Once reached, the heat-cycle light will cycle on and off with the heaters maintaining set temperature. Typically, the oven will need to cycle at a set temperature for a minimum of 20 minutes before it will achieve equilibrium and become temperature stable.

NOTE: The temperatures printed on the "quick-set" dial are designed to help quickly set a temperature to within a close proximity of the indicated dial temperature. Small rotational adjustments to the dial will likely be required to set a more precise temperature setting as measured against a reading from the door thermometer or other temperature measuring device.

Also, any degree of offset observed for a given temperature setting may be different for other temperature settings on the dial. The control dial is calibrated at the factory in the middle of the model's temperature range and is therefore most accurate in these middle-range temperatures. Over time, continuous use at a single temperature setting may require periodic re-adjustment as the contacts wear or as ambient temperatures change seasonally, or from air conditioning or heating. See also: "Temperature Dial Calibration".

Unit Safety Reset- If the unit exceeds its maximum temperature, an internal safety "snap disc" will cut power to the unit until it is manually reset. To do this, unplug the unit, lay it on its back and remove the bottom cover. Locate the round "snap disc" and push in the button to reset.

⚠ WARNING: RISK OF ELECTRICAL SHOCK. DISCONNECT UNIT FROM POWER SOURCE BEFORE REMOVING COVER. FAILURE TO COMPLY COULD RESULT IN SERIOUS INJURY OR DEATH.

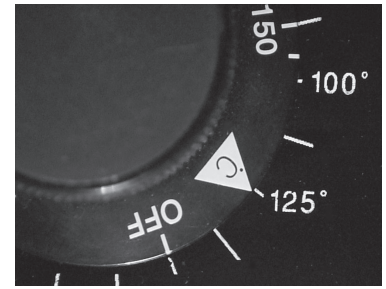


FIG. 4

Chamber Loading

Article or media processing times and temperature uniformity are largely dependent on load density and positioning. Load the oven so that the air circulation within the oven is not impaired. Here are some general guidelines:

Leave a space between multiple articles on a shelf as shown in (FIG 5). Avoid placing articles or media against or in close proximity to walls, allowing unrestricted air flow around articles and contributing to even and consistent heating.

Use of large solid trays, or foil on shelves severely limits the oven's ability to distribute heat evenly and uniformly. See (FIG 6). Since not enough heat rises within the chamber, door thermometer readings give false indication that temperature setting is too low. Higher temperature adjustments made as a result of these readings could overheat the lower-placed articles or media.

For best processing performance for a single item, adjust a shelf so that the item is centered in the oven chamber.

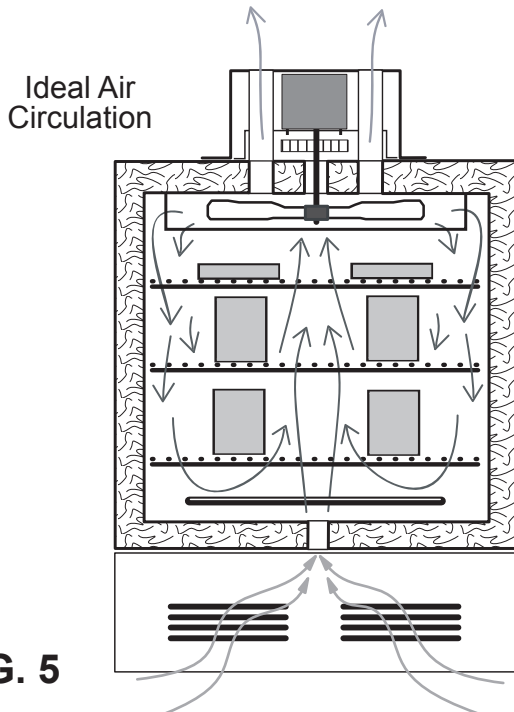


FIG. 5

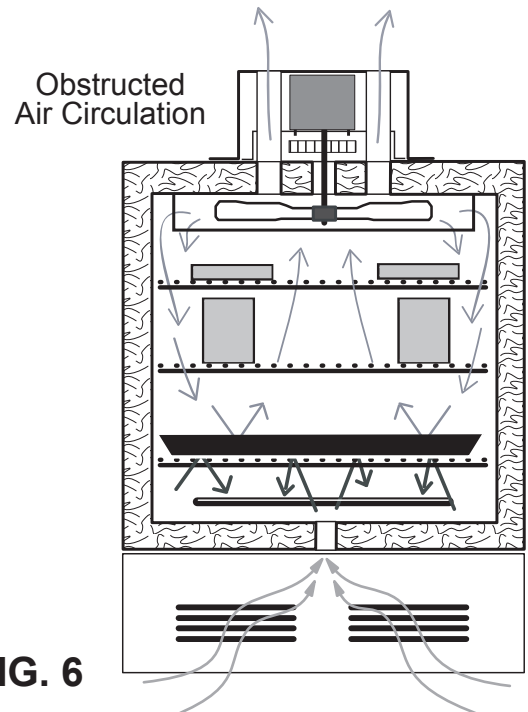


FIG. 6

Chamber Loading (continued)

Don't overload the unit with large (in quantity OR size), or high-density loads. This will restrict air flow and result in non-uniform processing and long or impossible "heat-through" times. To help determine a large load's suitability, use the set point recovery time (the time it takes for the temperature to recover to the original set temperature once load is placed) as a guide. To reduce recovery time, reduce load proportionally.

Care should also be taken to avoid placing items on the top-most shelf too close to the holes in the upper fan plenum (See upper plenum **FIG 1**). This will restrict the flow of air passing through the plenum and reduce the amount of pre-heated air being circulated through the entire chamber, causing uneven chamber temperatures.

NOTE: The set point of the thermostat or chamber temperature stability can be affected by changes in ambient (room) temperature and/or equipment running on the same electrical circuit. Take time to see how unit location or changes in room temperature from seasonal heating or air conditioning may influence the oven's set temperature. Isolate the oven on a separate circuit when possible.

The unit's minimum operating temperature is largely determined by ambient temperature. The unit can operate 25 degrees F (15 degrees C) above room temperature but temperature stability will be degraded. Stability improves appreciably for settings that exceed ambient by 40 degrees F (20 degrees C) or better.


Periodic Oven Maintenance


The AF series Lab Ovens are designed to be virtually maintenance free. But operational safety requires periodic cleaning and chamber temperature accuracy verification. Periodically check the rear air intake vents for dirt or dust build-up. Keep the intake & exit ports clear of obstruction, dust and dirt. Check the actual oven chamber temperature yearly against a known accurate temperature measurement device. Maintain temperature dial accuracy to within +/-15 degrees F of the control setting. Calibrate the control as necessary. To clean exterior and interior surfaces, use a damp cloth or an all-purpose cleaner. Avoid commercial oven cleaners.

Temperature Dial Calibration

Calibrating the actual chamber temperature to the temperature printed on the knob/dial, should only be done if the chamber temperature, as measured by a known accurate device, is off by more than +/- 15°F (8°C) from any indicated dial setting. The most accurate calibration possible for any analog-type control is at a single temperature setting. Therefore, calibrate the control at the most frequently used temperature.

To calibrate: Set dial temperature to a frequently used temperature then remove knob (allen screw). Be careful not to move the position of the shaft when removing the knob. Allow oven to achieve a stable temperature. Remove compression pin (newer models only) to access the trim-screw recessed in shaft center. Using a small blade driver, turn the trim-screw clockwise (CW) to adjust temperature down and (CCW) to adjust the temperature upward. Make only small 1/16th adjustments at a time. Allow the temperature to settle before making any subsequent adjustments. **NOTE:** Use pliers to keep outer shaft from turning when making trim-screw adjustments.

 Do not adjust control to operate above the unit's stated maximum operating temperature.

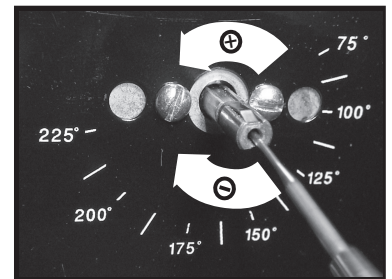

Technical Support

If you have any questions or need technical assistance, please contact Quincy Lab Tech Support at

Email: information@quincylab.com
Voice: 800-482-HEAT (4328)
Fax: 773-622-2282

Quincy Lab, Inc.
1925 N Leamington Ave
Chicago, Illinois 60639

Visit us on the web at www.quincylab.com

Limited Warranty



Quincy Lab, Inc. warrants to the original purchaser that this product will be free from defects in material and workmanship under normal use throughout the warranty period. The standard warranty period for this instrument is 18 months from date of shipment. The instrument warranty is supplemented by a three year warranty on the heating elements. Please refer to your invoice / shipping documents to determine the active warranty period. Warranty covers parts & labor (labor at factory only) and shipping costs for replacement parts.