




17120 Tye Street SE, Suite D * Monroe, WA 98272
Ph (360) 453-2090 Ph (800) 289-6753 Fx (360) 453-2091
Email: sales@awscientific.com www.wine-testing-supplies.com

TABLET TEST COLOR COMPARISON CHART FOR SUGAR IN WINE

- 1) Pipet $\frac{1}{2}$ ml of sample into test tube. (or use 10 drops from dropper supplied with some kits)
- 2) Drop a sugar test tablet (Clinitest) into the tube.
- 3) Watch the tube as a boiling reaction proceeds. If the color goes past orange to brown, the sample is too sweet to test and must be diluted before retesting.
- 4) Wait 15 seconds after boiling stops, then compare the color in the tube with this color chart. (Do not use Clinitest chart.)
- 5) Estimate the % sugar directly from the color chart.

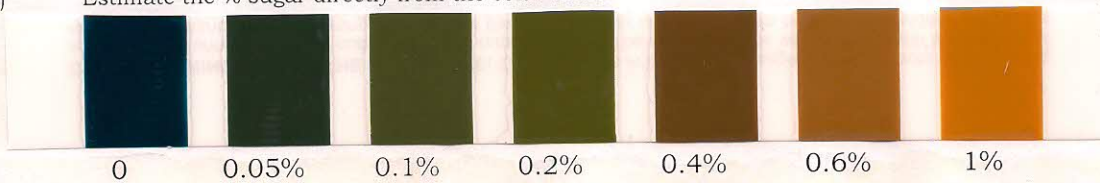

% RESIDUAL SUGAR PRESENT

(accuracy approx. +/- 0.05 to 0.1%)