

# Pilot-Scale Universal Soft Gel Capsule Machine

SGL207NextGen®









## Pilot-Scale Universal Soft Gel Machine

## SGL207NextGen®

The **SGL207NextGen®** Pilot-Scale Universal Soft Gel Machine has an operational speed of up to 5 RPM, 4 medicine plungers, and electronic gearing using servo motors. Its robust construction meets cGMP requirements. PLC and touchscreen provide more reliable control over production and process parameters for improved accuracy. Its design assists with easy cleaning, maintenance, assembly, and convenient changeover. It can be integrated with dryers, polishers, chilling systems, etc. Suitable for R&D, laboratories, nutraceutical industries, cannabis, and other facilities. **Available for both gelatin and veggie-based films** with our TES VGen closed boxes and our TES VGen metering system.

### **Features**

- Pilot-scale soft gelatin machine.
- Motion Control System: the most important mechanical gearing replaced with electronic gearing in synchronized loops, using servo motors.
- Die Roll dimensions: approximately 4" x 3".
- Operation speed 0.5 5 RPM.
- Control of all production and process parameters using Allen Bradley<sup>®</sup> programmable Logic Controllers (PLC) with touchscreen.
- Electrical panel and PLC incorporated within the machine.
- Sealed construction to prevent any possible rust and allows easy cleaning and maintenance.
- The outside covers of the machine are entirely made from stainless steel, designed according to the GMP/FDA requirements.
- CFR 21 Part 11 compliance software.

## **Mechanism**

## **Medicine Pump System**

- The medicine pump is independently driven by a servo motor and electronically synchronized with the die rolls.
- During a capsule size change, the operator inputs the number of die roll cavities using the touchscreen, and the medicine pump timing is done electronically (no gear change is required).
- Precise injection timing adjustments are done electronically using the touchscreen; The timing can also be adjusted during operation without stopping the machine. Since the timing can also be adjusted with such high accuracy, this improves on seal integrity of the capsule, which will increase capsule quality.
- The slide lubrication system for the pump is integrated into the drive mechanism.
- Four plunger standard pump.

#### **Spreader Box**

- Adaptable gelatin ribbon width that allows a minimum loss of ribbon net.
- Bronze construction.
- Gelatin level control is done electronically for gravity feeding from individual attached tanks; ground-level gelatin feeding is also available using a pressurized gelatin tank (optional).

#### **Casting Drums**

- "In-line" construction allows a straight gelatin ribbon transfer to the oil rolls and die rolls without any twisting.
- Water/Air-cooled cast drums.

## Pilot-Scale Universal Soft Gel Machine

## SGL207NextGen®

#### **Ribbon Lubrication & Drive System**

- Porous material for lubrication roller.
- Electronic metering pump and a precise distributor valve allow accurate ribbon lubrication.
- Easy and accurate oil quantity adjustment.
- Independent electrical drives allow fine adjustments of the ribbon tension on each side.
- Easy access from the operator's side to feed the ribbon on the oil rollers.

## **Mechanism**

#### **Die Roll Pressure System**

• Manually operated pneumatic valves for high and low pressure.

#### The Divider (Chute Assembly)

- Variable speed independent drive.
- Sealed construction.

### **Mangle Rolls**

• Electronic speed adjustment allows fine-tuning of the ribbon tension and helps the cutting process.

### **Wedge Lifting System**

• Air-actuated lifting with a safe middle position of the injection wedge.

#### **Machine Lubrication**

• Closed lubrication system with a dedicated pump for the Die Roll housing.

### **Soft Capsule Transfer Conveyor**

- Variable speed drive motor.
- Bi-directional.


## Set of aluminum die rolls (4" x 3") and change parts

- Specially coated injection wedge.
- Storage case.
- Die size per customer requirements.

# Pilot-Scale Universal Soft Gel Machine

# SGL207NextGen®

# **Specifications**

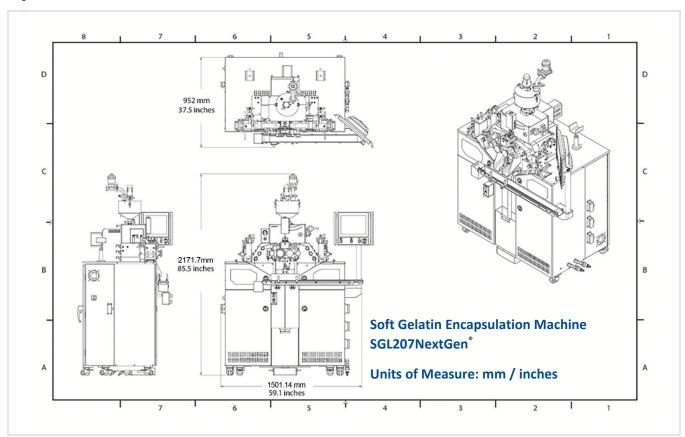

| Model No.                |   |
|--------------------------|---|
| SGL207NextGen®           | Soft Gel Capsule Machine Next Gen - 5RPM                                      |
| Specifications           |   |
| Dimensions (mm/inches)   | 1501.14mmL x 952mmW x 2171.7mmH (mm) / 59.1" L x 37.5" W x 85.5" H x (inches) |
| Weight (kg/lb)           | 900 kg / 1984 lbs.  |
| Machine Electrical Power | 15 KVA  |
| Compressed Air           | 100 PSI   |
| Chilled Water            | 3.8 GPM   |



**Technophar Global Headquarters**