

ⓘ When using this homepage, we recommend Chrome browser. Chrome can translate web

pages.

ABOUT US

PRODUCT

APPLICATIONS

FAQ

CONTACT

HIRANUMA APPLICATION DATA		Karl Fischer Titrator	Data No.	KF19	Feb. 10, 2021
Water contents	Sugars – KF Volumetry, granulated sugar Factor standardization with water standard				

1. Abstract

Water contents of granulated sugar could be determined by Karl Fischer volumetric titrator. In volumetric titration, titrant have a factor which means the amount to react with water per 1 mL of titrant. The factor is pre-determined before sample measurement and water content of sample is calculated from the consumed titrant volume by sample measurement.

This application introduces an example for the water determination in granulated sugar. Due to the low water content of this sample, a relatively large amount of sample must be added, but the solubility in methanol as titration solvent is not sufficient. Sugars have tendency to dissolve in formamide. Therefore a mixed solvent of formamide and methanol was chosen as the titration solvent. To make it easier to dissolve the sample, the titration cell was heated to 45 °C using an outer chamber for flowing warm water.

Since the water content of the granulated sugar is often a few hundred ppm, the titrant with low factor value, as 1 mg/mL was used. When pure water is used as the standard material of factor standardization for low factor titrant, the amount of pure water added should be as small as 10 to 20 mg, which makes accurate addition and weighing difficult. Therefore, water standard material which is the solution containing 1 % of water sealed in glass ampoule is used for the factor standardization.

2. Apparatus and Reagents

(1) Apparatus

Titration cell	: Titration cell with cooling chamber P/N D327510-1
	* This cell could not fit to the stirrer of AQV-300.
Powder funnel	: less than 12 mm outer diameter of leg
Syringe	: Glass, 10 mL

(2) Reagents

KF Titrant	: HYDRANAL Composite 1 (Honeywell)
methanol	: HYDRANAL Methanol (Honeywell)
Formamide	: HYDRANAL Formamide dry (Honeywell)
Titration solvent	: Mixed solvent of methanol and formamide at a volume ratio of 1: 2
Standard material	: Water standard 10 (certificate of analysis : 10.00 mg/g)

* Equivalent product can be substituted these reagents.

(3) Other

Heating device	: Temperature-controlled bath with circulation function
	* Commercial item

3. Procedure

3.1. Factor standardization

ⓘ When using this homepage, we recommend Chrome browser. Chrome can translate web

pages.

controlled bath, to heat the titration solvent.

[ABOUT US](#)

[PRODUCT](#)

[APPLICATIONS](#)

[FAQ](#)

[CONTACT](#)

(4) Open an ampoule of water standard 10 and collect the contents in a syringe. To prevent injury when opening the ampoule, handle it according to the instructions of standard material.

To make easier to collect, it is helpful to use packing box of standard material to stand the ampoule as shown by an example in Fig.3.2.

(5) Put the syringe on the balance. Record its read (S1 [g]). Inject sample from rubber septum of titration cell as shown in Fig.3.3.

(6) Start titration. Measurement parameter is shown in Table 4.1.

(7) Weigh the syringe again and record its read (S2 [g]). The difference of (S1-S2 [g]) is set as sample size.

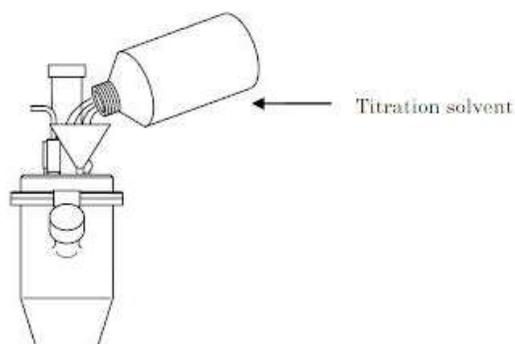


Fig.3.1 Preparation of the reagents

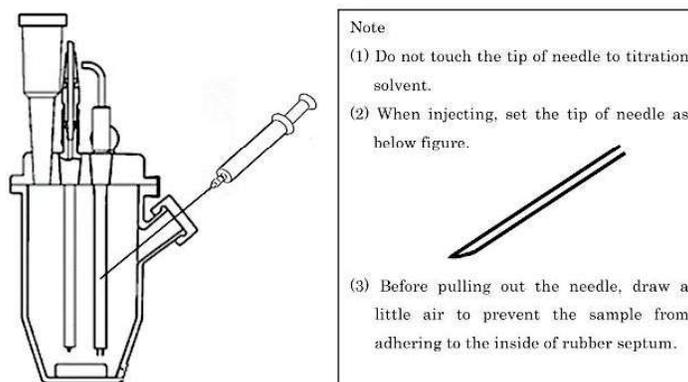


Fig.3.2 Example of water standard ampoule and its packing box.

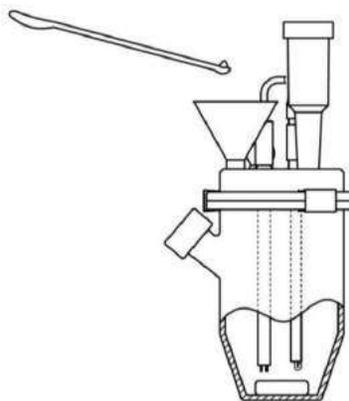


Fig.3.3. Injection of sample.

3.2. Sample measurement

(1) Fill 50 mL of titration solvent into the titration cell as shown in Fig.3.1.

ⓘ When using this homepage, we recommend Chrome browser. Chrome can translate webitrolled bath, to heat the titration solvent.

pages.

[ABOUT US](#)

[PRODUCT](#)

[APPLICATIONS](#)

[FAQ](#)

[CONTACT](#)

(5) Open the glass stopper of titration cell and to introduce the sample with powder funnel as shown in Fig.3.4.

(6) Start titration. Measurement parameter is shown in Table 4.2.

(7) Weigh the sample container, powder funnel and spatula again and record its read (S2 [g]). The difference of (S1-S2 [g]) is set as sample size.

* If the sample adheres to the spatula or funnel, include these in the tare of balance.

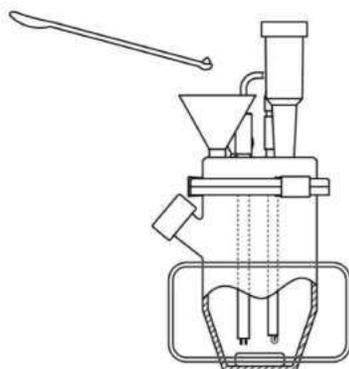


Fig.3.4 Introduction of sample with powder funnel

4. Parameters and results

Table 4.1 Parameters of factor standardization

Item			
Cal Mode	8:KF Factor(By Standard)		
Interval Time	30 sec		
Max Volume	20 mL	Sample Size Input	Every Time
Min Feed Vol.	0.01 mL	Blank Value	0 mL
S.Timer	0 min	E.P Detection	mV
KF Buret No.	1	Solvent	FM
STD H ₂ O Conc.	1.000 %	C.P Level	150 mV
KF Speed(OUT)	24 mL/min	E.P Level	4
KF Speed(IN)	24 mL/min	Auto Interval	0 g
Back Ground	ON	Auto Input	OFF

Table 4. 2 Parameters of sample measurement

Item			
Cal Mode	0:Sample weight (net)		
Interval Time	30 sec		
Max Volume	20 mL	Sample Size Input	Every Time
Min Feed Vol.	0.01 mL	Blank Value	0 mL
S.Timer	10 min	Unit Mode	AUTO
KF Factor	1.0853 mg/mL	E.P Detection	mV

ⓘ When using this homepage, we recommend Chrome browser. Chrome can translate web

ABOUT US	PRODUCT	APPLICATIONS	FAQ	FM
Back Ground	ON	Auto Interval		150 mV
				CONTACT
				4
				0 g

Table 4.3 Results of factor standardization

Sample	Meas. No.	Sample size (g)	BG (mL)	Titrant volume (mL)	Factor (mg/mL)	Statistics result	
	1	0.8749	0.020	8.13	1.0788	Avg.	1.0853 mg/mL
Water	2	1.0943	0.036	10.11	1.0863	SD	0.0051 mg/mL
standard	3	1.0697	0.052	9.87	1.0895	RSD	0.47 %
10 mg/g	4	1.0814	0.054	9.97	1.0906		
	5	1.0263	0.050	9.54	1.0815		

Table 4.4 Results of water content measurement in granulated sugar

Sample	Meas. No.	Sample size (g)	BG (mL)	Titrant volume (mL)	Water (mg)	Water content (ppm)	Statistics result	
	1	3.1130	0.096	0.81	0.775	249.0	Avg.	239.2 ppm
	2	3.1342	0.130	0.81	0.738	235.5	SD	9.2 ppm
Granulated	3	3.1568	0.119	0.82	0.761	241.1	RSD	3.86 %
sugar	4	6.2303	0.121	1.48	1.475	236.7		
	5	6.2351	0.132	1.56	1.550	248.6		
	6	6.2127	0.136	1.42	1.394	224.4		

5. Note

- (1) The blanking might become unstable when time has passed since the solvents of methanol and formamide were mixed. The cause could be due to the ammonia generated by mixing these solvents. Adding 3 g of benzoic acid to 50 mL of the titration solvent improves the unstable state of blanking.
- (2) In the measurement of granulated sugar, the dissolution time of the sample was set to 10 minutes in the parameter "S-Timer". If the sample becomes difficult to dissolve due to repeated measurements, replace it with a new solvent. In this report, the titration solvent was replaced with a new one every 3 sample measurements.
- (3) In the measurement of this report, the measurement condition "Background" is set to ON for background correction. The background is the consumed volume of KF titrant to keep the titrant solvent anhydrous during blanking. With this function, the amount of titrant consumed by the background is subtracted from the measurement result.

The reasons for using the background correction in this report are as follows:

In the measurement of this report, the experimental conditions are such that the blanking is relatively less stable than in the normal measurement, and about 0.01 mL/min of KF titrant is consumed in blanking. This is due to the following experimental conditions.

- (i) Low factor of KF titrant
- (ii) Formamide-containing titration solvent

ⓘ When using this homepage, we recommend Chrome browser. Chrome can translate web

pages.

trant consumed to maintain the anhydrous

[ABOUT US](#)

[PRODUCT](#)

[APPLICATIONS](#)

[FAQ](#)

[CONTACT](#)

and the background has a large effect on the measurement results, so the correction is required.

le measurement is approximately 1 mL,

Keywords: Karl Fischer, Volumetric titration, Granulated sugar, Formamide, Heating



[ABOUT US](#)

[PRODUCT](#)

[APPLICATIONS](#)

[FAQ](#)

[CONTACT](#)

[TERMS OF USE](#)

[PRIVACY POLICY](#)

