

Toll Free: 866-702-0260 Phone: 203-433-4329 Fax: 203-208-0664

www.phoenixsongsbio.com

FL-HCC Spheroid Dissociation Medium

Catalog Number: 35301-030

For Dissociation of

Human Fibrolamellar Hepatocellular Carcinoma (FL-HCC) Spheroids

Compete Serum-Free FL-HCC Spheroid Dissociation Medium is formulated for dissociation of human FL-HCC spheroids. FL-HCC Spheroid Dissociation medium enzymes necessary for the dissociation of the spheroids with optimal survival of the cells and is filtered to ensure sterility, frozen and maintained in -80°C ultralow freezer until ready for use. The medium is shipped on dry ice and when thawed, the medium is complete and ready to use for dissociation of human FL-HCC spheroids.

Handling FL-HCC Spheroid Dissociation Medium

- 1. FL-HCC Spheroid Dissociation medium (Cat.# 35301-030) is complete and ready to use:
 - a. Store in -80°C ultralow freezer until ready for use
 - For optimal performance place frozen bottle of media in 4°C refrigerator to thaw overnight. The next day it will be ready to use.
 - c. You may also thaw in 37°C waterbath
 - i. Place frozen bottle of media in waterbath
 - ii. Invert bottle every 5-10 minutes until only a small ice ball remains then remove from the waterbath
 - d. DO NOT leave media in 37°C waterbath for longer than 5 minutes as this may result in degradation of the growth factors, proteins and glutamine in the media
 - e. Store thawed media in 4°C refrigerator in the dark
 - f. Use FL-HCC Spheroid Dissociation medium within one month
 - g. DO NOT re-freeze FL-HCC Spheroid Dissociaiton Medium

Reference:

Model of Fibrolamellar hepatocellular carcinomas reveals striking enrichment in cancer stem cells. 2015. Tsumekazu Oikawana, Eliane Wauthier, Timothy A. Dinh, Sara R. Selitsky, Andrea Reyna-Neyra, Guido Carpino, Ronald Levine, Vincnzo Cardinale, David Klimstra, Eugenio Gaudio, Domenico Alvaro, Nance Carrasco, Praveen Sethupathy and Lola M. Reid. Nature Comm. 6, Article 8070. DOI: 10.1038

Use Restrictions: The Products and components of the Products are to be used for RESEARCH PURPOSES only and are not for use in humans or animals. No rights expressed or implied are conveyed by the sale of the Products to make, have made, offer to sell or sell the Products alone or in combination with any other products or to use for commercial purposes without a sub-license from PhoenixSongs.