

HOW TO LOOK AT STOMATA

WITH A MICROSCOPE!

YOU WILL NEED:



Leaf



Clear Nail Polish



Tape



Microscope



1. APPLY NAIL POLISH!

Put a thin layer of clear nail polish on the underside of the leaf.



2. LET IT DRY

Wait 10-15 minutes until the polish is fully dry.



3. PEEL IT OFF

Carefully peel off the dried nail polish from the leaf.



4. STICK ON TAPE

Place the polish peel on a piece of clear tape.



5. LOOK CLOSELY!

Put it under the microscope and observe the stomata!



WHAT YOU'LL SEE:

STOMATA UNDER THE MICROSCOPE

