

NOCOBOX



■ ■ Manufactured in France by OxyPharm according to ISO 13485.

The **Nocobox** are culture dishes with nutrients to highlight germs on surfaces. They are used to verify the result of disinfection from an airborne surface disinfection system. They contain a nutrient that promotes the growth of bacteria and therefore highlights their presence.

STORAGE

Store the boxes in a cool (below 30°C) and dry place. Store out of sunlight. Length of validity in the original closed packaging: 4 months.

Before



After



IMPORTANT

Elimination: collection and incineration by specialised companies.

INSTRUCTIONS FOR USE

1st sampling before disinfection

- 1 Open the box carefully (do not touch the test area).
- 2 Apply 10 seconds of light pressure on the surface to be tested.
- 3 Close the box.
- 4 Leave it at development temperature (20°C - 25°C) for 48 hours.
- 5 Read the results.

2nd sampling after disinfection

- 6 Repeat steps 1 to 5 after disinfection using new boxes.
- 7 Compare and observe the effectiveness of the treatment.

REFERENCES

PACKAGING

REFERENCE

Boxes of 10

5000.020