

Instruction for Use

*alpha*Clean Proteinase K

Molecular biology grade

For general laboratory use.
For *in vitro* use only.

REF

831007



50 mg



1 Package Contents

Molecular biology grade Proteinase K, 1 x 50 mg lyophilized.

Proteinase K is an endolytic protease that cleaves peptide bonds at the carboxylic sides of aliphatic, aromatic or hydrophobic amino acids. Proteinase K is a broad-spectrum serine protease that is active under denaturing conditions and stable at high temperatures. The smallest peptide to be hydrolyzed by this enzyme is a tetrapeptide.

2 Product Specifications

Source:	<i>Tritirachium album</i>
Activity:	≥2.5 U/mg (Casein, 37°C)
RNase Activity:	Not detectable
DNase Activity:	Not detectable
Inhibitors:	DFP, Hg ²⁺ , PMSF

One unit is defined as the amount of enzyme that will liberate 1 µmol of tyrosine per minute at 37°C, pH 7.5.

3 Transport and Storage

The *alpha*Clean Proteinase K must be stored at +18 to +25°C. If properly stored, the enzyme is stable until the date of expiry printed on the label.

Reconstituted Proteinase K solution has to be aliquoted. Aliquots stored at ≤ -18°C are stable for 12 months.

4 Procedure

When using with *alpha*Clean Pure RNA/DNA Kit or *alpha*Clean Mag RNA/DNA Kit reconstitute Proteinase K in 2.5 ml PCR grade water (final concentration: 20 mg/ml).

5 Abbreviations and Symbols



Catalog number




Batch code



Contains sufficient for <n> test



Content

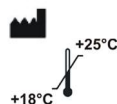

Manufacturer



Use by YYYY-MM

Limit of temperature



Consult instructions for use



gerbion gmbH & Co. KG

Remsstr. 1
70806 Kornwestheim
Germany
phone: +49 7154 806 20 0
fax: +49 7154 806 20 29
e-mail: info@gerbion.com

Distributor

Mikrogen GmbH

Floriansbogen 2-4
D-82061 Neuried
Germany

phone: +49 (0) 89-54801-0

fax: +49 (0) 89-54801-100

e-mail: mikrogen@mikrogen.de

www.mikrogen.de